



Sterile, alcohol based ready to use solution for surface disinfection in cleanrooms

perform® sterile alcohol IPA

At a glance:

- Fast, effective and residue free
- Sterile (gamma irradiated) and microbial filtrated (0.2 µm)
- Ready to use in a handy, sterile 500ml spray system
- Optional with WFI (water for injections)
- Colour and perfume free
- Efficacy tested to European norms

Applications

Disinfection of surfaces and equipment / devices in cleanrooms, in particular for class A and B cleanroom areas, according to:

- Annex 1 of EU Guide to good manufacturing practice: Manufacture of sterile medicinal products, March 2009
- FDA Guidance for Industry: Sterile drug products produced by aseptic processing – current good manufacturing practice; U.S. Department of Health and Human Services, Food and Drug Administration; September 2004.

Instructions for use

Apply perform® sterile alcohol IPA undiluted to objects and surfaces and allow it to work, ensuring that all areas are moistened completely. Remove any excess solution with a disposable cloth. Do not use to treat surfaces that are particularly sensitive to alcohol, such as acrylic glass. perform® sterile alcohol IPA is also available with WFI.

Microbiological effectiveness

perform® sterile alcohol IPA is: bactericidal • levurocidal • fungicidal • virucidal (BVDV, Vaccinia).

Contact times

bactericidal	1 Min.
levurocidal	1 Min.
fungicidal	15 Min.
virucidal (BVDV, Vaccinia)	30 Sec.



Tested according to EN 1276, 1650, 13697

Product data

Composition: 100 g contain: 63.1 g Propan-2-ol (equates 70% (v/v))

Chemical / physical data:

Appearance:	colourless liquid
Density at 20 °C:	0.88 g/cm ³
Flash point (DIN 51755):	22 °C

Labelling according to EC Directives

Xi:	Irritant
R10:	Flammable
R36:	Irritating to eyes
R67:	Vapours may cause drowsiness and dizziness
S23:	Do not breathe vapour
S24/25:	Avoid contact with skin and eyes
S26:	In case of contact with eyes, rinse immediately with plenty of water and seek medical advice
S51:	Use only in well ventilated areas

perform[®] sterile alcohol IPA

Special notes

Use disinfectants safely. Always read the label and product information before use.

The sterile 500 ml spray system offers the following advantages:

- Convenient size for storage.
- Ergonomically shaped spray pump for ease of use.
- Remains sterile after opening.

The bottles are sterile (with irradiation indicator) and double-wrapped individually.

perform[®] sterile alcohol IPA comprises alcohol of pharmaceutical quality in accordance with the Ph. Eur.

With correct handling and storage the in-use product life corresponds to the use by date printed on the product label.

Wearing of protective gloves (e.g. butyl rubber) is recommended.

Forms of delivery/packaging units

	Container sizes	Packaging units	Art. No.
Standard	500 ml spray bottle	10 x 500 ml	136406
WFI	500 ml spray bottle	10 x 500 ml	121101

Environmental information

schülke manufactures products economically and with advanced, safe and environmentally friendly production processes while at the same time maintaining our high quality standards.

Expert opinions and information

Please visit our website for an overview of all available literature/ reports on perform[®] sterile alcohol IPA: www.schuelke.com

For further information please contact your local representative or local offices.



Schülke & Mayr GmbH is certified according to DIN EN ISO 9001, DIN EN ISO 14001 and DIN EN ISO 13485 (Reg.-No. 004567-MP23) and has a validated environmental system in accordance with the Eco Audit Regulation (Reg.-No. DE-150-00003).

Schülke & Mayr GmbH
22840 Norderstedt, Germany
Phone +49 (0) 40 - 521 00 - 0
Fax +49 (0) 40 - 521 00 - 247
www.schuelke.com
mail@schuelke.com

Schülke & Mayr UK Ltd.
1 Jenkin Road
GB-Sheffield S9 1AT
Phone +44 (0) 1142 - 54 35 - 00
Fax +44 (0) 1142 - 54 35 - 01
mail.uk@schuelke.com

Schülke & Mayr Ges.m.b.H
Seidengasse 9
1070 Wien, Austria
Phone +43 (0) 1 - 523 25 01 - 0
Fax +43 (0) 1 - 523 25 01 - 60
office.austria@schuelke.com