



Sterile, ready to use 70 % (v/v) IPA impregnated wipes for surface disinfection in cleanrooms.

## perform® sterile wipes IPA

### Our Plus

- fast, effective and residue free
- pouch of 20 ready to use alcohol impregnated wipes
- sterile (X-Ray irradiated) and double wrapped
- suitable for cleanrooms grade A/B, ISO class 5/6
- low particle release
- colour and perfume free
- efficacy tested according to European norms (EN)

### Application areas

Disinfection of surfaces and equipment / devices in cleanrooms, in particular for class A and B cleanroom areas, according to:

- Annex 1 of EU Guide to Good Manufacturing Practice Manufacture of sterile medicinal products, March 2009.
- FDA Guidance for Industry: Sterile drug products produced by aseptic processing - current Good Manufacturing Practice; U.S. Department of Health and Human Services, Food and Drug Administration, September 2004.

### Instructions for use

Apply perform® sterile wipes IPA to objects and surfaces and allow to take effect, ensuring that all areas are wiped completely. The product dries off quickly. Do not use to treat surfaces that are particularly sensitive to alcohol, such as acrylic glass. The wiping reach depends on the properties of the surface to be disinfected and the ambient conditions.

### Microbiological efficacy

Efficacy	Concentration	Contact time
<b>bactericidal</b> <i>EN1276, EN13697</i>	ready-to-use	1 min.

Efficacy	Concentration	Contact time
<b>yeastcidal</b> <i>EN1650, EN13697</i>	ready-to-use	1 min.
<b>virucidal against enveloped viruses</b> <i>in accordance with DVV (German Registered Association for Combating Viral Diseases)//RKI Guideline</i>	ready-to-use	30 sec.



# perform<sup>®</sup> sterile wipes IPA

## Product data

### Composition:

100 g solution contains the following active substances: 63.1 g Propan-2-ol (equates 70% (v/v)).

### Chemical-physical data

Color	colourless
Density	ca. 0,88 g/cm <sup>3</sup> / 20 °C
Flash point	22 °C / Method : DIN 51755 Part 1
Form	liquid
pH	Not applicable
Viscosity, dynamic	> 2,5 mPa*s / 20 °C / Raw material

## Special advice

**Use disinfectants safely. Always read the label and product information before use.**

The alcohol used within the perform<sup>®</sup> sterile wipes IPA comprises alcohol of pharmaceutical quality in accordance with the Ph. Eur.

EU-0023656-0002 1-2

## Information for order

Item	Delivery form	Item no.
perform <sup>®</sup> sterile wipes IPA 20 pc bag	20/Carton	on request

These products are not available in every country. For more information please contact our local subsidiary or distributor.

## Related Products

- perform<sup>®</sup> sterile alcohol EP
- perform<sup>®</sup> sterile alcohol IPA
- perform<sup>®</sup> sterile alcohol IPA (WFI)
- perform<sup>®</sup> sterile dry wipes
- perform<sup>®</sup> sterile mix dry wipes
- perform<sup>®</sup> sterile premium dry wipes

## Environmental information

schülke manufactures products economically and with advanced, safe and environmentally friendly production processes while at the same time maintaining our high quality standards.

## Expert opinion and information

Please visit our website for an overview of all available literature/reports on the product: [www.schuelke.com](http://www.schuelke.com). For individual questions:

Customer Sales Service

Phone: +49 40 52100-666

E-Mail: [info@schuelke.com](mailto:info@schuelke.com)



Schülke & Mayr GmbH holds a Manufacturer's Authorisation according to sect 13 para 1 German Drug Law and Certificates of GMP Compliance for medicinal products.

**schülke Headquarters**  
Schülke & Mayr GmbH  
Robert-Koch-Str. 2  
22851 Norderstedt, Germany  
Phone +49 40 - 52100 - 0  
Fax +49 40 - 52100 - 318  
[www.schuelke.com](http://www.schuelke.com)  
[info@schuelke.com](mailto:info@schuelke.com)