



Sterile, ready to use 70 % (v/v) IPA impregnated wipes for surface disinfection in cleanrooms.

perform® sterile wipes IPA

Our Plus

- fast, effective and residue free
- pouch of 20 ready to use alcohol impregnated wipes
- sterile and double wrapped
- suitable for cleanrooms grade A/B, ISO class 5/6
- low particle release
- colour and perfume free
- efficacy tested according to European norms (EN)

Application areas

Disinfection of surfaces and equipment / devices in cleanrooms, in particular for class A and B cleanroom areas, according to:

- Annex 1 of EU Guide to Good Manufacturing Practice Manufacture of sterile medicinal products, March 2009.
- FDA Guidance for Industry: Sterile drug products produced by aseptic processing - current Good Manufacturing Practice; U.S. Department of Health and Human Services, Food and Drug Administration, September 2004.

Instructions for use

Apply perform® sterile wipes IPA to objects and surfaces and allow to take effect, ensuring that all areas are wiped completely. The product dries off quickly. Do not use to treat surfaces that are particularly sensitive to alcohol, such as acrylic glass. The wiping reach depends on the properties of the surface to be disinfected and the ambient conditions.

Microbiological efficacy

Efficacy	Concentration	Contact time
bactericidal <i>EN1276, EN13697</i>	ready-to-use	1 min.

Efficacy	Concentration	Contact time
yeastcidal <i>EN1650, EN13697</i>	ready-to-use	1 min.
virucidal against enveloped viruses <i>in accordance with DVV (German Registered Association for Combating Viral Diseases)//RKI Guideline</i>	ready-to-use	30 sec.



perform[®] sterile wipes IPA

Product data

Composition:

100 g solution contains the following active substances: 63.1 g Propan-2-ol (equates 70% (v/v)).

Special advice

Use biocides safely. Always read the label and product information before use.

For all of the related safety information please refer to the Safety Data Sheet. To request the Safety Data Sheet please E-mail: tech.uk@schuelke.com

The alcohol used within the perform[®] sterile wipes IPA comprises alcohol of pharmaceutical quality in accordance with the Ph. Eur.

Information for order

Item	Delivery form	Item no.
perform sterile wipes IPA 20 ST BT	20/Carton	121103

Environmental information

schülke manufactures products economically and with advanced, safe and environmentally friendly production processes while at the same time maintaining out high quality standards.

Expert opinion and information

Please visit our website for an overview of all available literature/reports on all schulke products online at <http://www.schuelke.com/>

For any queries you might have about the product or using it please call 0114 2543500 or E-mail mail.uk@schuelke.com



Schülke & Mayr GmbH holds a Manufacturer's Authorisation according to sect 13 para 1 German Drug Law and Certificates of GMP Compliance for medicinal products.

schülke UK
Schülke & Mayr UK Ltd
Cygnet House, Meadowhall
Sheffield, S9 1AT
Phone +44-114-254 35 00
Fax +44-114-254 35 01
mail.uk@schuelke.com

schülke Headquarters
Schülke & Mayr GmbH
Robert-Koch-Str. 2
22851 Norderstedt, Germany
Phone +49 40 - 52100 - 0
Fax +49 40 - 52100 - 318
www.schuelke.com
info@schuelke.com